

Packing List for 2pcs/ 8pcs cartridge kit

	DNA cartridge Kit		RNA cartridge Kit		Volume (μl)	
	2pcs	8pcs	2pcs	8pcs	2pcs	8pcs
Alignment Marker	●	●			100*2	100*8
Alignment Marker 5x Lower Mark			●	●	100*2	100*8
Separation Buffer	●	●			50	250
10x Separation Buffer			●	●	15	50
Dilution Buffer	●	●			15	50
10x Dilution Buffer			●	●	8	25
Mineral Oil	●	●	●	●	8	25
Accessory Package: Buffer Tray, PCR Vial and Dropper	●	●	●	●	-	-



BiOptic



Website: www.biopic.com.tw

Email: info@biopic.com.tw

Tel: +886-2-2218-8726

Fax: +886-2-2218-8727

Address: (23141) 4F, No. 108-3, Minquan Rd., Xindian Dist., New Taipei City, Taiwan (R.O.C.)

Unpacking Guide



1. Open gel cartridge calm shell container and lift the cover to the bottom.



2. Take cartridge out vertically and remove the stickers from the top.



3. Take out the pin and push it straight into the hole of the cartridge cap.



**Cartridge is ready to use.
Enjoy it !!!**

5. Remove the pin and clean the cartridge tip by tissue.



4. Push the pin down to the end then move up and down for 2-3 times.

