

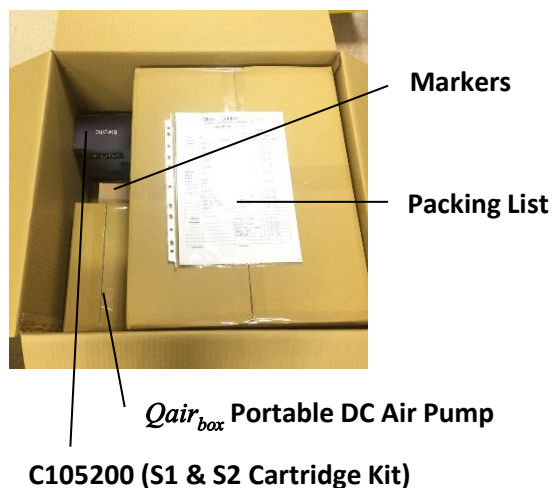


# Getting Started with *Qsep*<sub>100</sub>

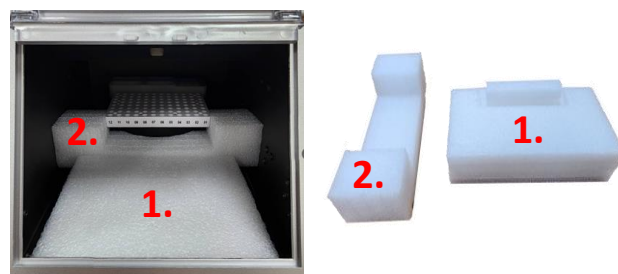

## Installation Steps

1. Unpack the *Qair*<sub>box</sub> and securely tighten the air tube with the connector.
2. Plug the power cord into the *Qair*<sub>box</sub>. Ensure the label on the power cord matches *Qair*<sub>box</sub>.
3. Remove the protective foams **before** powering on the *Qsep*<sub>100</sub>.
  - Retain the protective foams for securing the tray holder during transportation
4. Plug the power cord and USB cord into the *Qsep*<sub>100</sub>. Ensure the label on the power cord matches *Qsep*<sub>100</sub>.

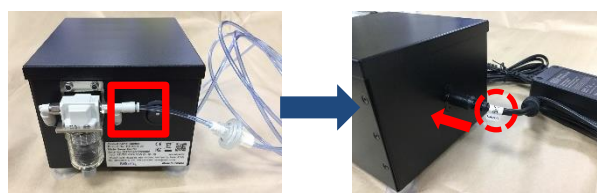

5. Connect the USB cord to the operation computer.




6. Plug the other end of the air tube into the *Qsep*<sub>100</sub> and turn on both the *Qair*<sub>box</sub> and *Qsep*<sub>100</sub>.



## Packing Procedure

1. Remove buffer tray, samples, and cartridge.
2. Click "Lock" to initiate the movement of the tray holder to its initial position.

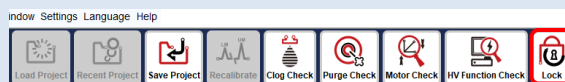

3. Turn off the *Qsep*<sub>100</sub> and push the tray holder to the very back.



4. Place the protective foams to secure the tray holder.



- Insert the foam into the gap to ensure stability.