

# Getting start with Qsep<sub>1</sub>-Plus



Packing List  
包裝清單

Markers  
標準品

C105200 (S1 & S2 Cartridge Kit)  
C105200 (S1 & S2 卡匣組)

Qsep<sub>1</sub>-Plus Power Adaptor  
Qsep<sub>1</sub>-Plus 電源供應線

Qsep<sub>1</sub>-Plus



Mini Centrifuge  
迷你離心機

SD Card & Software Key  
SD記憶卡及軟體金鑰

C105200 (S1 & S2 Cartridge Kit)  
C105200 (S1 & S2 卡匣組)

Qair<sub>box</sub> & Power Adaptor  
Qair<sub>box</sub> 及電源供應線

Ethernet Cable  
網路線

Sample Trays  
(8, 12 and 16+3 wells)  
樣品盤  
(8孔, 12孔和 16+3孔)

## Install Qsep<sub>1</sub>-Plus 安裝 Qsep<sub>1</sub>-Plus

1. Unpack Qair<sub>box</sub> and tighten the air tube with connector  
將 Qair<sub>box</sub> 開箱並把氣管插入轉接頭




Connector

2. Plug power cord into Qair<sub>box</sub>  
(\* Please check the Qair<sub>box</sub> label on the power cord)  
請先將 Qair<sub>box</sub> 接上電源  
(\* 請確認電源線上的標籤是屬於 Qair<sub>box</sub>)



Label

3. Remove the protecting foam before turning on Qsep<sub>1</sub>-Plus  
開啟 Qsep<sub>1</sub>-Plus 前務必移除保護泡棉

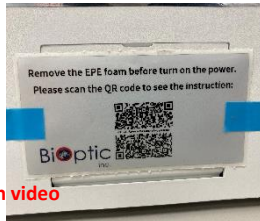


\* Scan the QR code for instruction video



4. Plug the power cord, Ethernet cable and the other side of the air tube into Qsep<sub>1</sub>-Plus and turn on the Qair<sub>box</sub> (Please make sure the label on the power cord is Qsep<sub>1</sub>)  
將電源供應線, 網路線及氣管線接上 Qsep<sub>1</sub>-Plus 並開啟 Qair<sub>box</sub>  
(請確認電源線上的標籤是屬於 Qsep<sub>1</sub>)



Your Qsep<sub>1</sub>-Plus is READY TO USE  
隨時可以使用您的 Qsep<sub>1</sub>-Plus

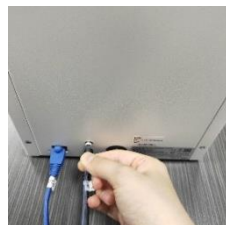

5. Connect the other side of the Ethernet cable to the operation computer.  
將網路線另一端連接您操作的電腦



**\*DO NOT switch on the instrument immediately after powering off  
Wait at least 5 seconds**